



HOW TO CONTROL CUSTOM PROFILE START IN ESI SOFTWARE

When you have a custom profile and want the ESI software to start the profile from the beginning and not from anywhere in the file, you may follow the following steps.


1. Add your OB1 to the ESI instruments list.

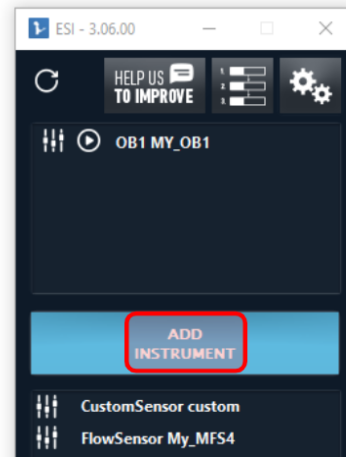


2. Load your custom profile

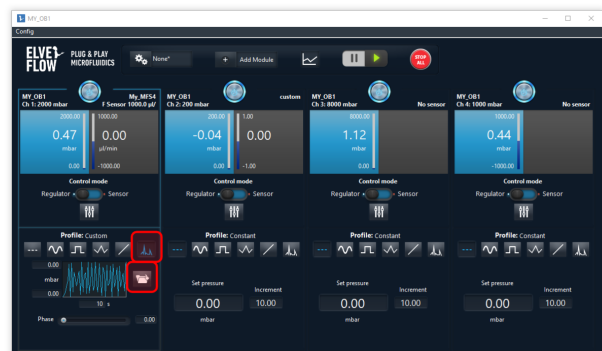


3. Save the Configuration

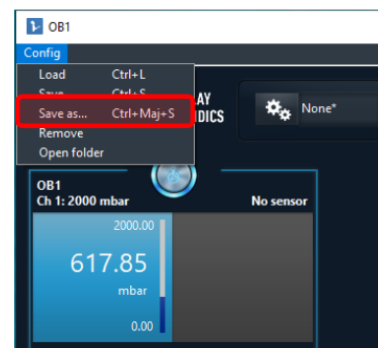

4. Open Sequencer

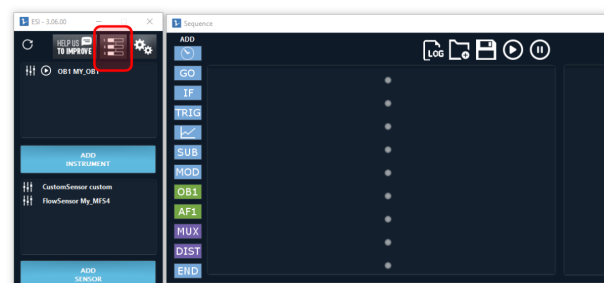

5. Create this sequence and save it (.sq file).

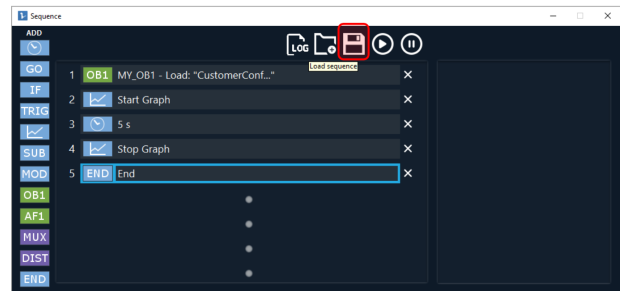


6. Open your sequence file (.sq) with Windows Notepad

7. Change the value of the “Unsynchronized” line from “False” to “True”, and save the file.



8. Load the sequence file on the sequencer

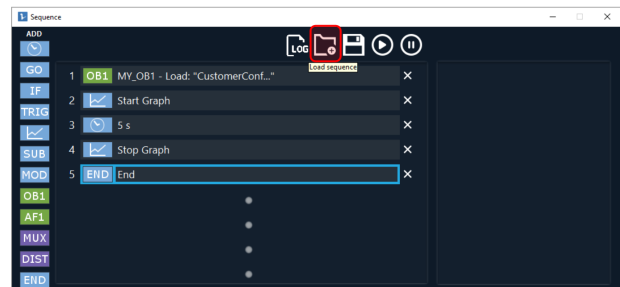

9. Run the program

